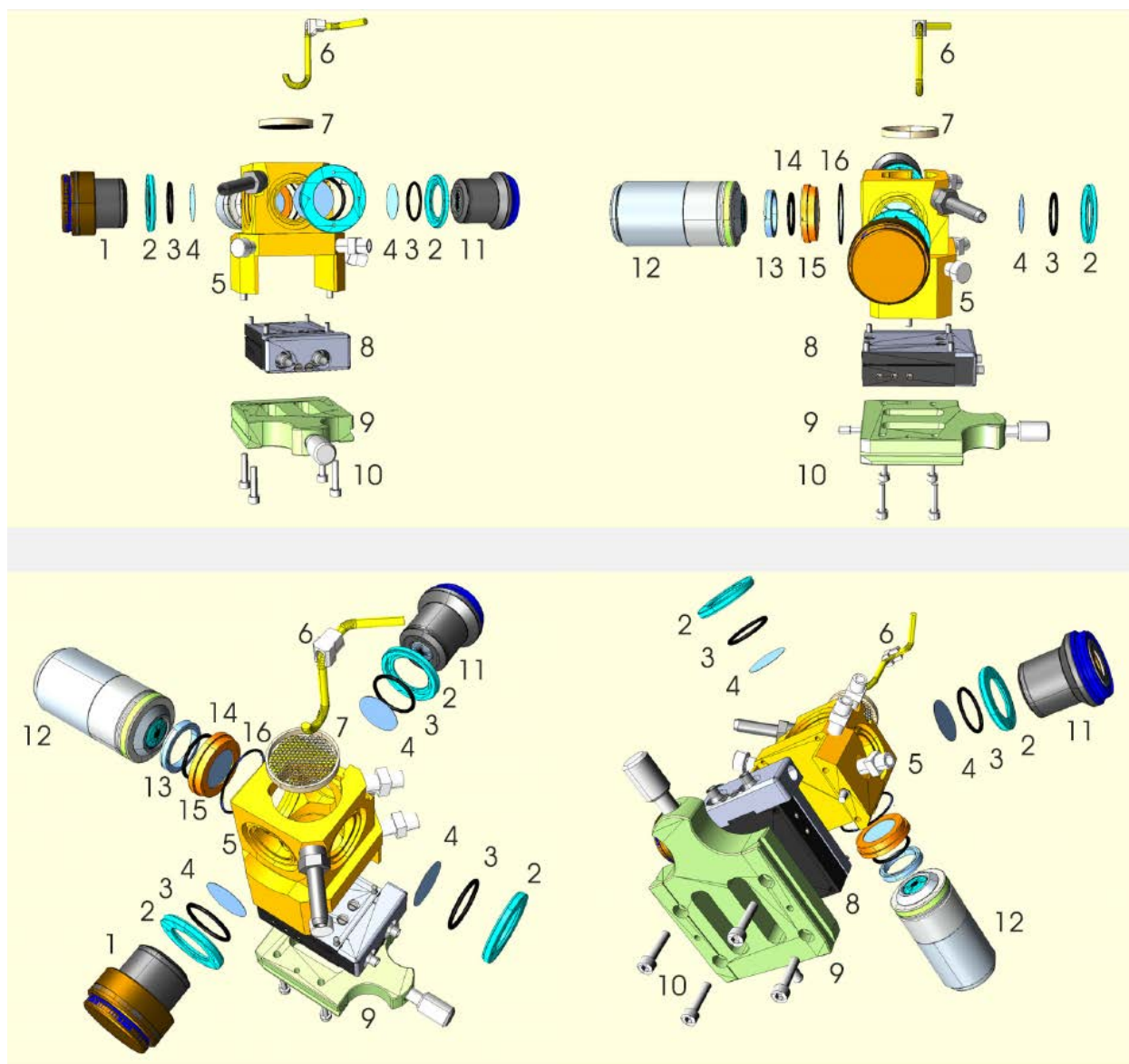
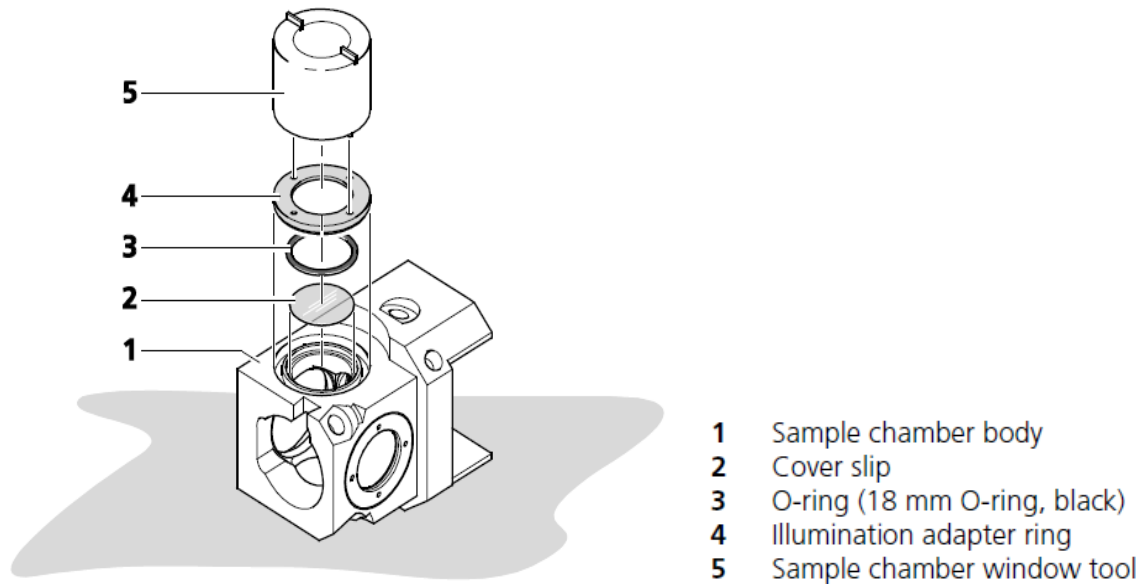
Sample Chamber 的安装

- | | | | | |
|----|---|----|--|--|
| 1 | Illumination objective, right-hand side, with cover ring (to be removed before use) | 12 | Detection objective | |
| 2 | Illumination adapter ring (screwed, for use with a cover glass) | 13 | <i>Air objective only:</i> Retaining ring (screwed, for use with objective 5x) | |
| 3 | Sealing ring (18 mm O ring, black rubber; order no. 0520-489) | 14 | <i>Air objective:</i> Sealing ring (18 mm O ring, black rubber; order no. 0520-489) | |
| 4 | Cover glass (order no. 0515-968) | 15 | <i>Water immersion objectives:</i> Sealing ring (15 mm O ring, black rubber; order no. 0515-968) | |
| 5 | Sample chamber, main body | 16 | <i>Air objectives:</i> Air retaining ring with cover glass (order no. cover glass 0515-968) | |
| 6 | PT100 temperature sensor | | <i>Water immersion objectives:</i> Water retaining ring without cover glass | |
| 7 | Strainer | | | |
| 8 | Heating block / Peltier element | | | |
| 9 | Base plate | | | |
| 10 | Fixing screws for carriage and heating block | | | |
| 11 | Illumination objective, left-hand side | | | |
| | | | 16 | Sealing ring (O ring, blue rubber, order no. 0520-488) |

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
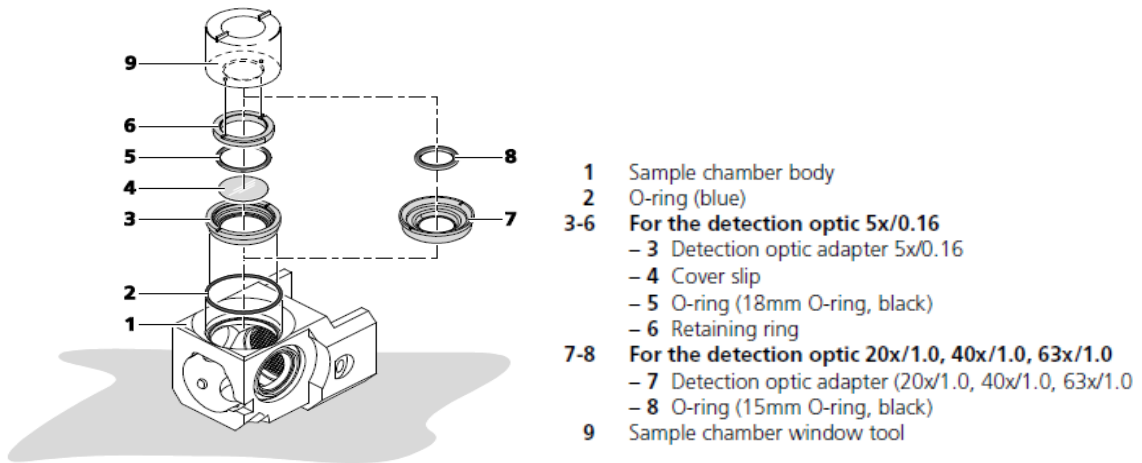
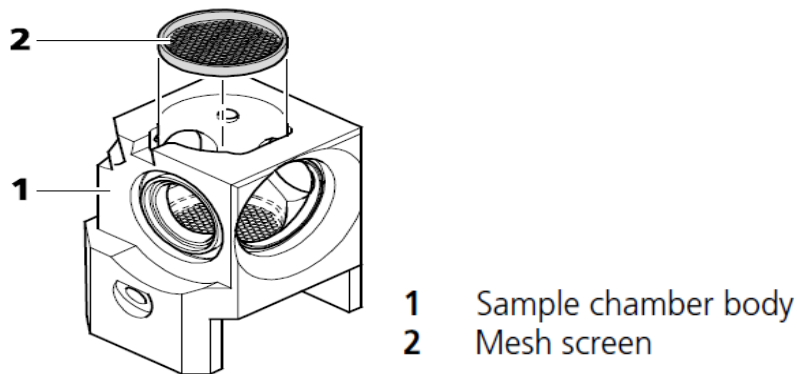
2. 安装 mesh screen 和 Sample Chamber Dove Tail Slide

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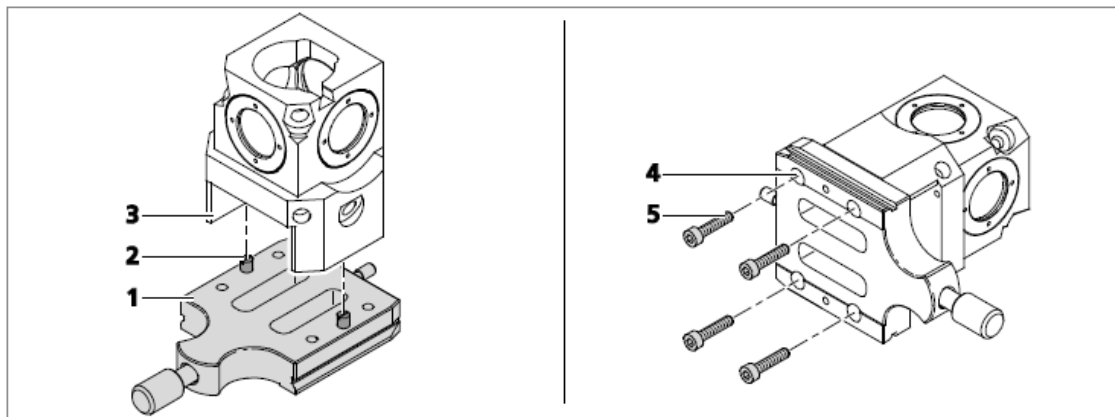
- 1 Sample chamber dove tail slide
- 2 Positioning pins
- 3 Sample chamber body
- 4 Screw-holes
- 5 Screws (diameter = M3, length = 14 mm)

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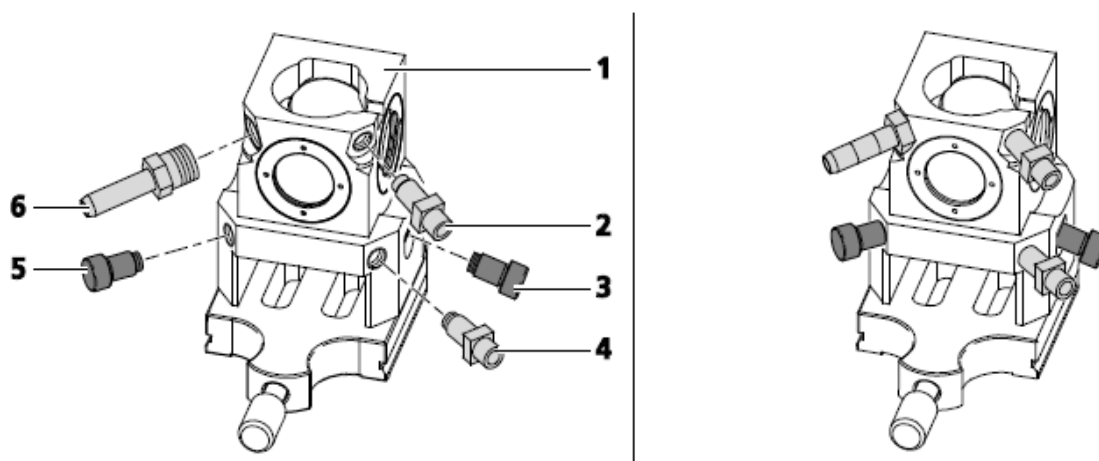3. 安装 Drain connector, Luer-Lock connector 和 Blind Plug

- 1 Sample chamber body with dove tail slide
- 2 Luer-Lock-connector (white); position recommend for CO₂ incubation
- 3 Blind plug (black)
- 4 Luer-Lock-connector (white); position recommend for syringe based perfusion
- 5 Blind plug (black)
- 6 Drain connector (grey)

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
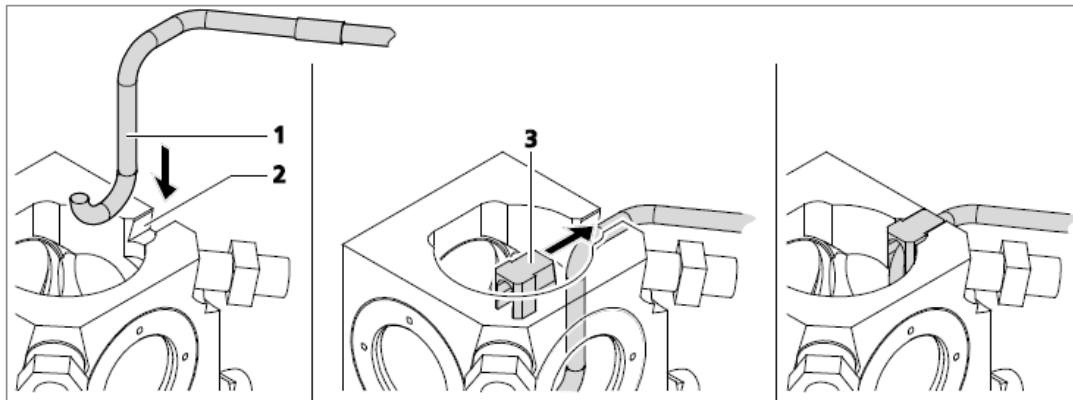
- 1 Temperature sensor
- 2 Sample chamber body groove
- 3 Temperature sensor sleeve

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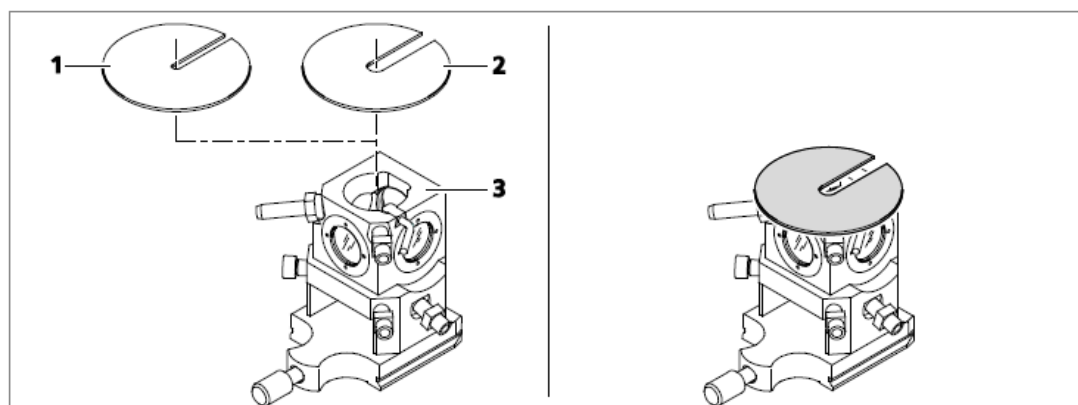- 1 Cover for capillaries (3 mm opening)
- 2 Cover for syringes (7 mm opening)
- 3 Sample chamber body with dove tail slide

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